



The final date for return of results is indicated on the accompanying request/report form

Contact details:

The Organisers - FEPTU
Public Health England
61 Colindale Avenue,
London, NW9 5EQ, UK.
Fax: +44 (0) 20 8200 8264
Tel: +44 (0) 20 8327 7119
e-mail: foodeqa@phe.gov.uk



Results should be returned on-line using the web-based reporting system. Go to: www.phe-ega.org.uk. You will need your log-in details for this process. Please contact us if you experience problems with the on-line system.

www.gov.uk/government/collections/external-quality-assessment-ega-and-proficiency-testing-pt-for-food-water-and-environmental-microbiology#public-health-scheme

Public Health Scheme - Instruction Sheet for food samples

The LENTICULE[®] discs are provided in screw-cap plastic vials (with a desiccant). Each LENTICULE disc rests on the filter support containing the desiccant. The **desiccant** should be orange in colour. Please contact us if this is not the case. Each LENTICULE disc requires reconstitution by a process of rehydration and dispersion prior to examination, as described below.

Refer also to the Safety Data Sheet: <https://www.gov.uk/government/publications/safety-data-sheet-for-lenticules>

Storage:

- Store the samples at **-20 ± 5°C** on receipt.
- Allow the LENTICULE discs to reach ambient temperature (5 - 10 minutes) **before** reconstituting in diluent.

Reconstitution:

- Open the sample container and transfer the LENTICULE disc into approximately **200mL of nutrient broth** that has been allowed to reach ambient temperature.
- Leave to stand at ambient temperature for 10 -12 minutes to rehydrate.
- Tighten the cap of the bottle and shake to disperse the micro-organisms.

Examination:

- Each reconstituted sample is equivalent to 200 gram of food.
- Undertake the sample examinations within 45 minutes of reconstitution.
- Examine in accordance with your routine procedures.

Return of Results:

- Return your results on-line using the web-based reporting system.
- The distribution will be closed on the deadline date for return of results.
- Results cannot be accepted after the distribution has closed.